

Introducing the Labplas SANI-STICK

- Ideal for surface sampling
- Allows to reach hard-to-reach areas
- Easier handling
- One time use
- Made to avoid sponge wedging
- Plastic sponge handle with a strong grip
- Comes with a 2.5 mil.in sterile sampling bag; 4.5 x 9 in
- **TWIRL'EM** bag with leak proof closure system
- Sterilized by gamma irradiation
- 100 units per case (10 packs of 10 units)
- **Choice of:**
 - Neutralizing agent
 - Neutralizing agent DE



Sani-stick user guide

STEP 1	STEP 2	STEP 3	STEP 4	STEP 5
Remove the tear-off strip and carefully take the sani-stick by the handle.	The wet sponge is already on the handle; swipe it on the determined surface.	Press on the bottom of the handle to gently release the sponge in the sterile sampling bag.	Twirl the bag three times and close it with the tabs.	Identify the sample on the write-on area with a permanent marker.

Code	Description	mil.in	micr.	in x in	mm x mm	oz	ml	Packaging
KSS-61310-NE	Sponge on handle with neutralizing buffer	2.5	63	4.5 x 9	114 x 229	15	450	Cs/100 (10 x 10)
KSS-61310-DE	Sponge on handle with DE neutralizing buffer	2.5	63	4.5 x 9	114 x 229	15	450	Cs/100 (10 x 10)

- STERILITY CERTIFICATE IN EVERY BOX/FDA GRADE PLASTIC
- 100% REPLACEMENT GUARANTEE ON ANY DEFECTS / CONTACT YOUR LABPLAS DISTRIBUTOR FOR FREE SAMPLES

CORPORATE HEADQUARTERS

1950 Bombardier Street
Ste-Julie, Quebec
Canada J3E 2J9

PHONE -CUSTOMER SERVICE

NORTH AMERICA
1.450.922.6946
OUTSIDE NORTH AMERICA
1.450.649.7343
Fax
1.450.649.3113
labplas@labplas.com

